

Herolab

Safety

The safety of our customers is very important to us. For this reason all our centrifuges have following safety features:

1. The stainless steel chamber has an additional inner guard ring.
2. Our APRI[®] system (Automatic Positive Rotor Identification) recognises the rotor itself and checks for the speed setting. In case the speed is selected too high for a particular rotor the machine will automatically reduce the speed setting to the maximum speed possible for the rotor in this centrifuge.
3. If any imbalance is detected an automatic cut-off will be activated.
4. There is a 4-point-lid fixing. The lid lock is motorised. There is also automatic lid opening and an emergency lid-lock release.
5. For all our centrifuges there is an overheat control.